

Understanding Small Molecule Biomarker Patterns by Targeted and Non-Targeted Metabolomics using LC/MS/MS

Henri Snijders

Applied Biosystems

Thessaloniki, 30 October 2008




Outline

- Non-targeted Metabolomics
- Targeted Metabolomics
- Novel Workflows for Lipidomics



Human Metabolic Pathways

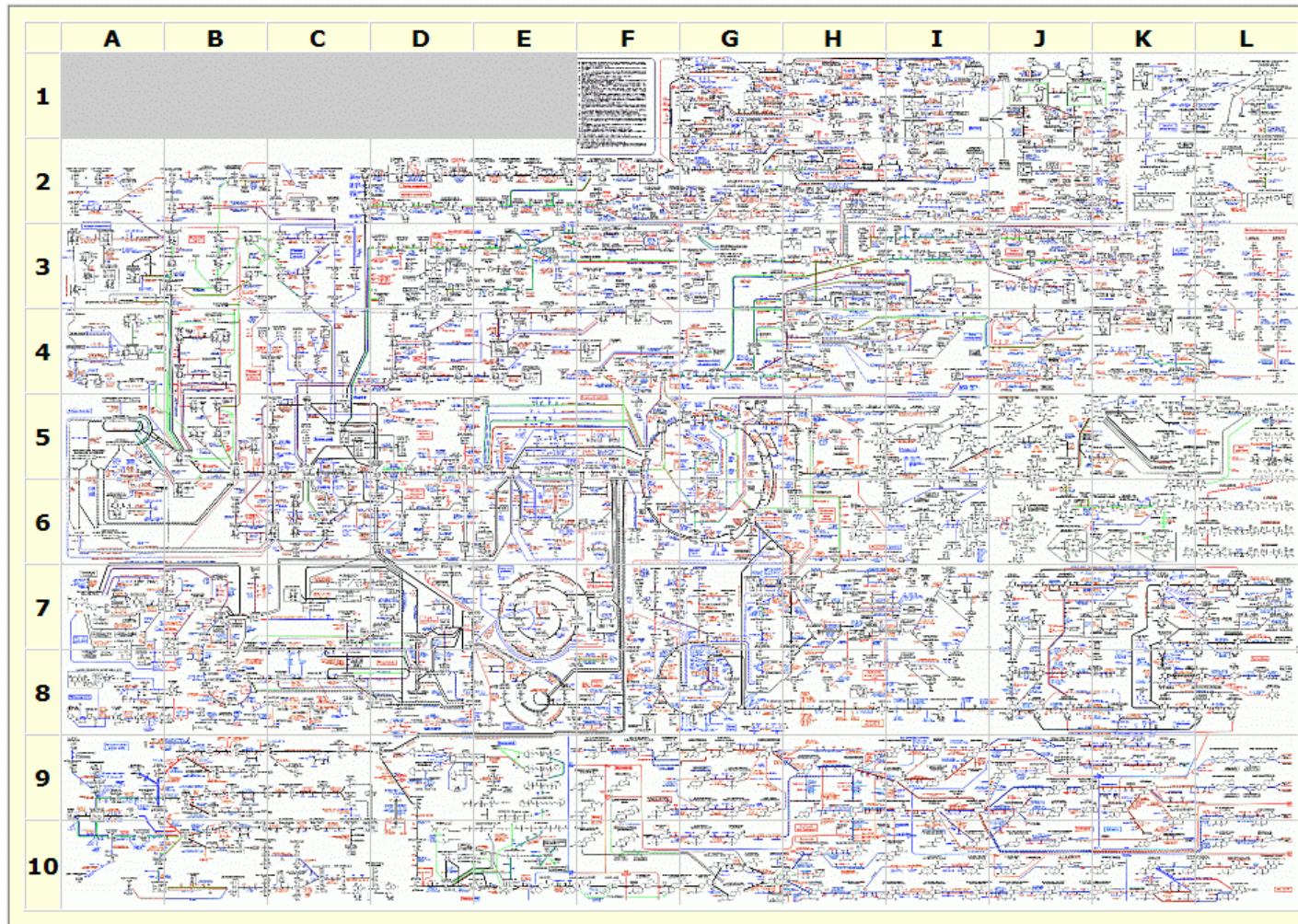



Metabolomics Workflows

| | | Application | | |
|----------|--------------|---|--------------|----------|
| | | Discovery | Pre-Clinical | Clinical |
| Workflow | Targeted | <p>Hypothesis-driven – have discovered important markers and need to monitor their presence and concentration</p> <ul style="list-style-type: none"> Acquire LC/MS data Monitor for known masses of interest Determine how these masses varied over multiple samples via stat analyses Find samples that are different (diseased, etc.) <p style="text-align: right; color: magenta;">4000 QTRAP/QTRAP 5500 QUAN</p> | | |
| | Non-Targeted | <p>Non-hypothesis-driven – not sure what you are looking for, screening large numbers of samples to look for patterns – diseased vs. non-diseased</p> <ul style="list-style-type: none"> Acquire LC/MS data <ul style="list-style-type: none"> – All unknowns or sample vs control (diseased vs nondiseased) Statistical analysis (PCA, HCA) to look for differences in samples Find samples that are different Identify what masses account for differences between samples Identify compounds <p style="text-align: right; color: purple;">QSTAR/ Qual</p> | | |



QSTAR[®] Elite LC MS/MS System

- Latest generation QqTOF platform
- Provides high resolution, accurate mass data (2 ppm and 15,000 resolution)
- 4 orders linear dynamic range in TOF mode
- Capable of acquiring 20 MS spectra/second and MS/MS on up to 7 precursor ions/second
 - Compatible with UPLC and other high throughput chromatography systems





Metabolomics Experiments

- What's really needed to for a metabolomics experiment?
 1. Must be able to find groupings in data (i.e. diseased samples must separate from non-diseased)
 2. Must be able to find potential biomarkers (unique masses that cause groups to separate)
 3. Must be able to determine elemental formula for potential biomarkers
 4. Must be able to identify potential biomarkers (assign name/structure to mass)
- What tools are best for this process?
 - What are the most important capabilities of hardware and software for a metabolomics experiment?
 - How important are scan speed, resolution, mass accuracy, isotope ratios, statistical analysis technique...?



Finding Groupings in the Data


- Typically acquire full-scan MS data and process using vendor-specific or general 3rd party statistical analysis tools



MarkerView™ 1.2 Software

Industry-leading Metabolomics Software

- Software for metabolomics and biomarker profiling experiments
- Finds and aligns peaks
- Performs statistical analysis to find potential biomarkers
 - T-Test
 - PCA
 - PCA-DA
- Can export data to any 3rd party stats package for additional processing
- View trends plots and link back to MS and MS/MS data
- Create IDA inclusions lists to acquire MS/MS data on selected compounds for further analysis
- Generate customized Word reports that can be shared amongst non-MarkerView users

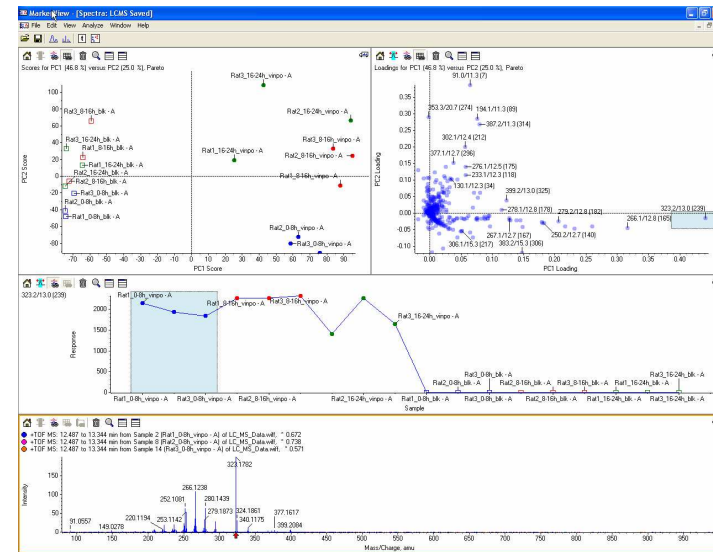



Finding Groupings in the Data

- Assume that if there are differences in the samples, some statistical package will enable you to find these differences
- However, can your data acquisition techniques affect whether you can see correct differences and groupings within data?
- Many metabolomics researchers moving towards high speed chromatography (UPLC) for faster separations and higher peak capacities
- What is the affect of these narrow LC peaks (1-2 seconds wide) on your data and your ability to correctly interpret data?
 - ... well, this depends on your MS scan speed



Finding Groupings in the Data

Importance of Scan Speed

- Can a slow-scanning instrument be used for metabolomics?
 - It depends....
- QSTAR[®] Elite System can acquire data at a rate of up to 20 spectra/second
- QTRAP[®] 5500 System has scan speed up to 20,000 amu/s and MRM dwell times down to 2 ms
- What happens to your ability to accurately interpret metabolomics data if you are acquiring only 1 spectrum/second and using high speed chromatography?




Experimental

- Samples:
 - Urine samples were collected in the morning from 3 different male and 3 different female subjects and frozen until analysis
 - A 1:10 dilution in mobile phase was done prior to analysis
 - Samples were divided into 9 aliquots, resulting in 9 replicate injections from each of 6 different subjects
- LC:
 - Chromatographic separation was performed using a PE 100 series autosampler and micro pumps
 - A gradient from 95% to 5% aqueous at 250 $\mu\text{L}/\text{min}$ through a BetaBasic C18 50 x 2.1 mm, 3 μm column was done over a 3 minute analysis
- MS:
 - A QSTAR® Elite System (Applied Biosystems|MDS Sciex) was used for collection of MS data over a 1000 amu mass range and all detection was performed in positive ion mode using TurbolonSpray® Source
 - Dynamic AutoCalibration was used to maintain mass accuracy throughout the data collection
 - Data acquisition was performed using several different scan speeds, ranging from 1 to 20 spectra/second
- Data Processing:
 - MarkerView™ Software



Typical TIC and XIC

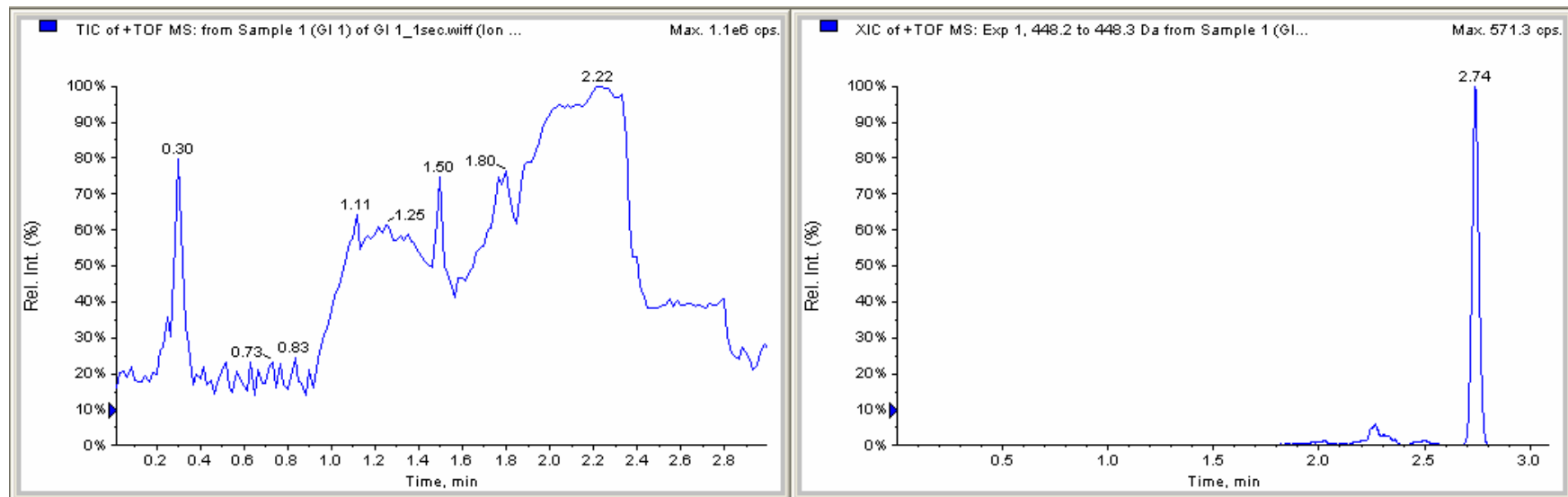


Data acquired at 20 spectra/second

XIC of m/z 448 showing the 4-second wide peaks typical from this experiment



Data Points Per Peak at different scan speeds

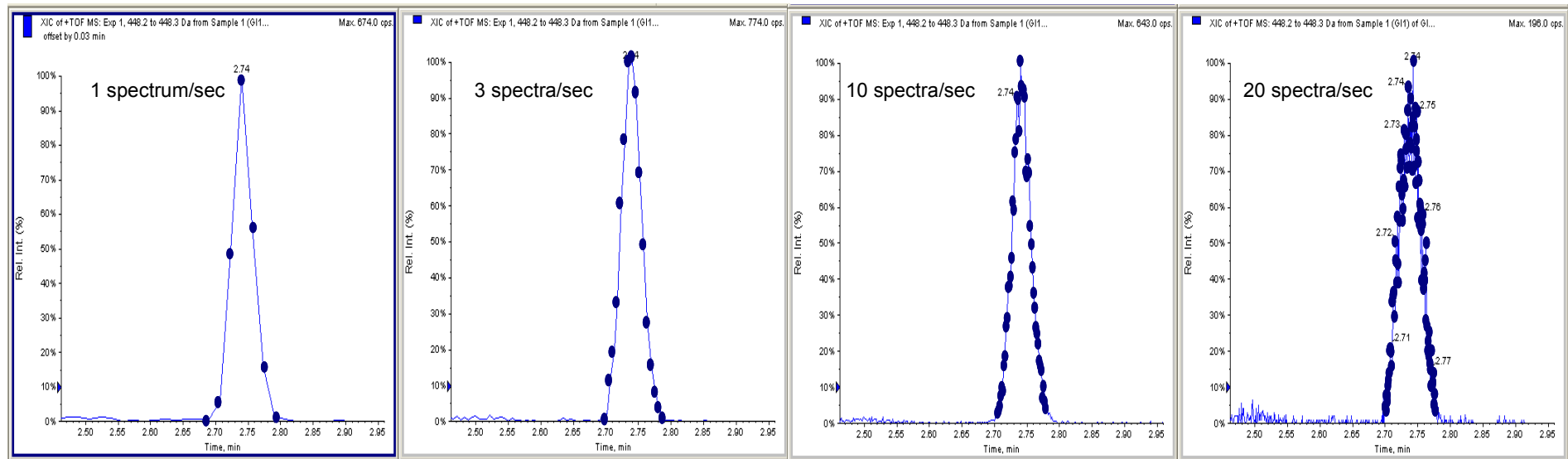

- At an acquisition rate of 1 spectrum/sec, only 3 data points above FWHH are acquired
- At an acquisition rate of 20 spectra/sec, ~40 MS spectra were acquired




PCA of data acquired at 1 vs 20 spectra/second

1 spectrum/second

20 spectra/second

- At 1 spectrum/second, there appear to be 3 distinct groups in the data
- At 20 spectra/second, there appear to be 6 distinct groups in the data
- Which is correct? Are there 6 groups, one for each individual, or 3 groups, one for male, one for female, and one for other?

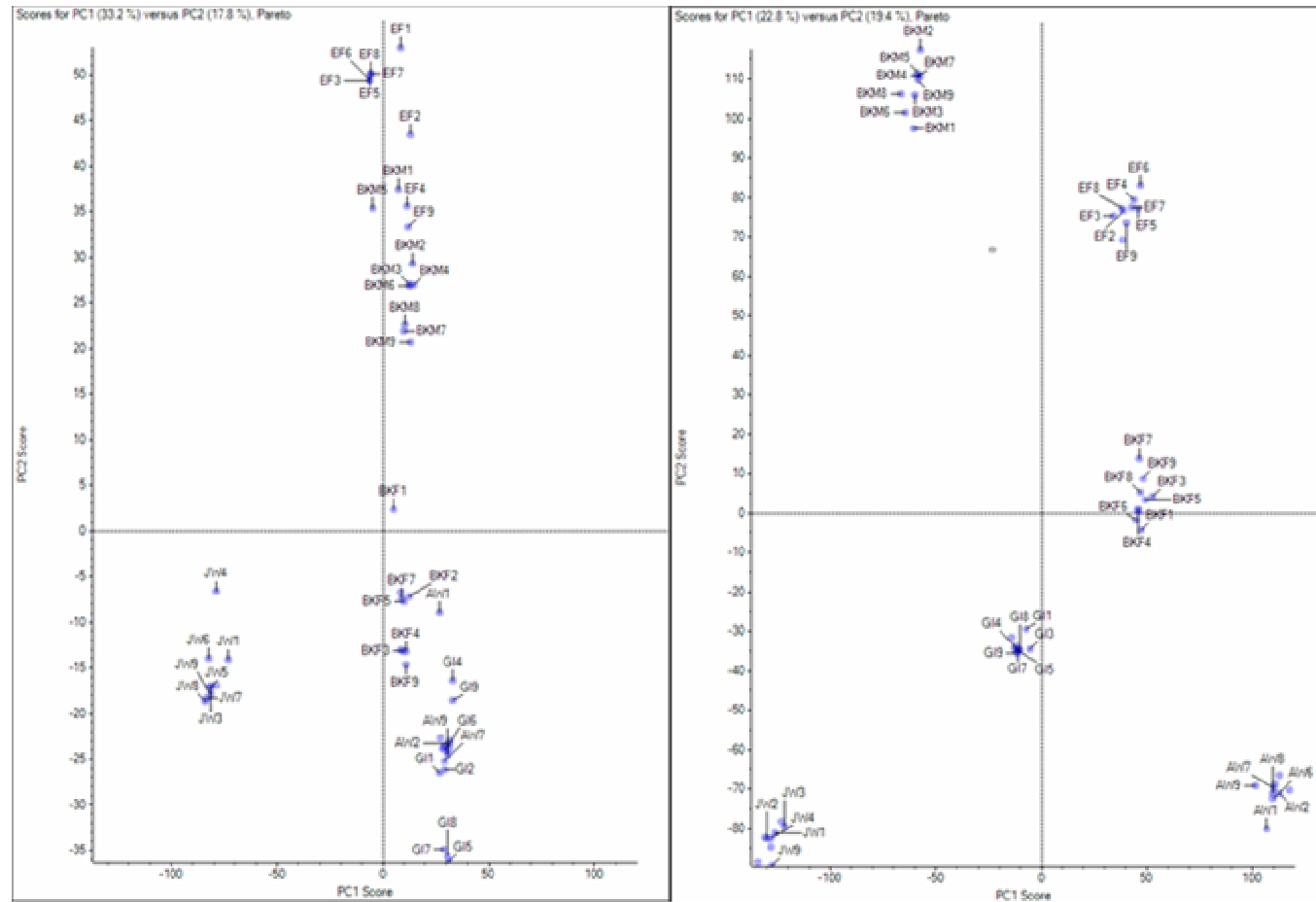




PCA of data acquired at 1 vs 20 spectra/second

1 spectrum/second

20 spectra/second

- Once colors are assigned to correct groups, becomes clear that the data acquired at 1 spectrum/second is very misleading
- Interestingly, at 1 spectrum/second, even the most basic distinction between men and women is incorrect
 - EF, BKF and JW are females
 - GI, AW, BKM are males

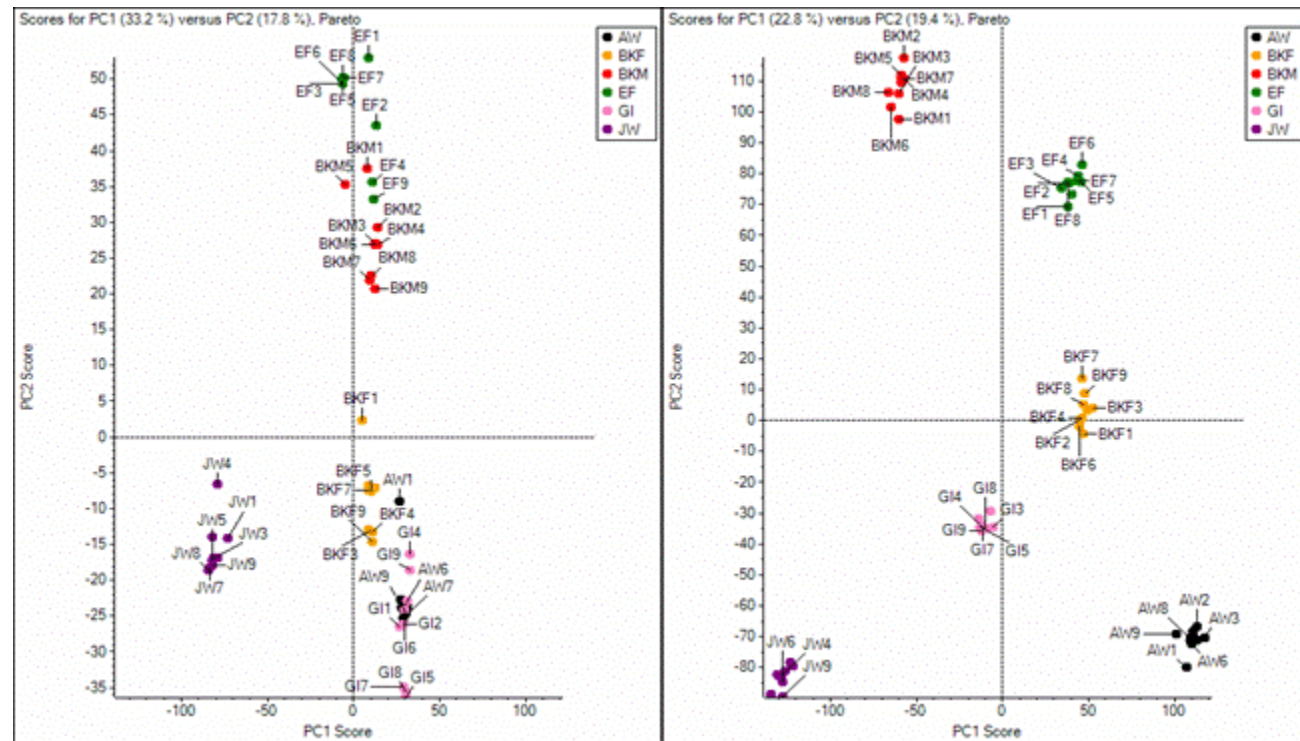

Data acquired at 1 spectrum/second is incorrect – shows only 3 groups within the data




Finding Potential Biomarkers

- Assume you have performed experiment correctly and have sufficient mass spec scan speed for the LC system being used and have reproducible data
- You are able to correctly group data and find m/z values that separate samples into groups



Determining Elemental Formula

- You now have a list of m/z values (along with retention times) for potential biomarkers
- Generally, to determine the structure of a compound, the first step is to determine the elemental formula
- Standard elemental calculators have the drawback that they often produce lists of 10's to even 1000's of possible formulae for a given molecular weight, even using a 5 ppm mass tolerance
- What's the solution?
 - Going to instruments with better mass accuracy precision?
 - Maybe not provided you can make more intelligent use of the data that you already have

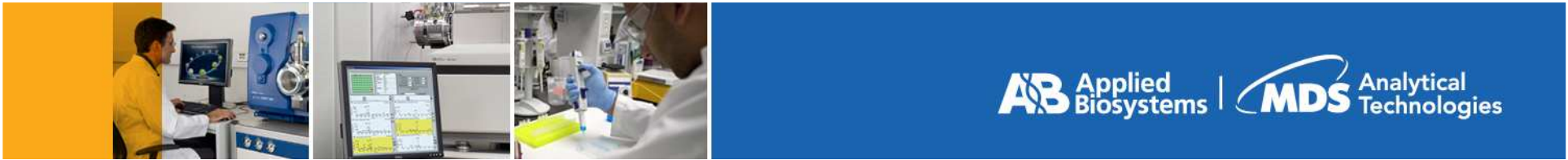

Formula Finder on the QSTAR® Elite


Simplifies compound identification

- Make use of all of the information available to make a better decision
- Uses a compound's molecular weight, isotopic pattern as well as applying 'chemical logic' to determine potential elemental formulae
 - Nitrogen rule (If the protonated pseudo molecular ion gives an odd mass then the number of nitrogen atoms is even)
 - Maximum number of double bond equivalents
 - Number of heteroatoms allowed
- Provides much shorter list of candidate formulae with the correct formula scoring much higher than standard elemental calculators




Reduction of Chemical "Space" using Chemical Sense

| Number | Name | Exact Mass | Formula |
|--------|--------------------|------------|---------------|
| | | MH+ | M |
| 1 | Detromethorphan | 272.2008 | C18H25NO |
| 2 | Quinoxifen | 308.0039 | C15H8Cl2FNO |
| 3 | Alprazolam | 309.0907 | C17H13ClN4 |
| 4 | Phenylbutazone | 309.1597 | C19H20N2O2 |
| 5 | Bestatin | 309.1808 | C16H24N2O4 |
| 6 | Benoxinate | 309.2172 | C17H28N2O3 |
| 7 | Lorazepam | 321.0192 | C15H10Cl2N2O2 |
| 8 | Omeprazole | 346.1219 | C17H19N3O3S |
| 9 | Tamoxifen | 372.2321 | C26H29NO |
| 10 | Haloperidole | 376.148 | C21H23ClFNO2 |
| 11 | Buspirone | 386.2556 | C21H31N5O2 |
| 12 | Clindamicin | 425.1877 | C18H33ClN2O5S |
| 13 | Loperamide | 477.2309 | C29H33ClN2O2 |
| 14 | Ketoconazole | 531.156 | C26H28Cl2N4O4 |
| 15 | Rescinnamine | 635.2969 | C35H42N2O9 |
| 16 | Bromocriptine | 654.2291 | C32H40BrN5O5 |
| 17 | Erythromycin - H2O | 716.4579 | C37H65NO12 |
| 18 | Roxithromycin | 837.5318 | C41H76N2O15 |
| 19 | Tylosin | 916.5264 | C46H77NO17 |
| 20 | Bromazepam | 316.00790 | C14H10BrN3O |
| 21 | Oxfendazole | 316.07500 | C15H13N3O3S |
| 22 | Oxycodone | 316.15430 | C18H21NO4 |
| 23 | Fendiline | 316.20590 | C23H25N |

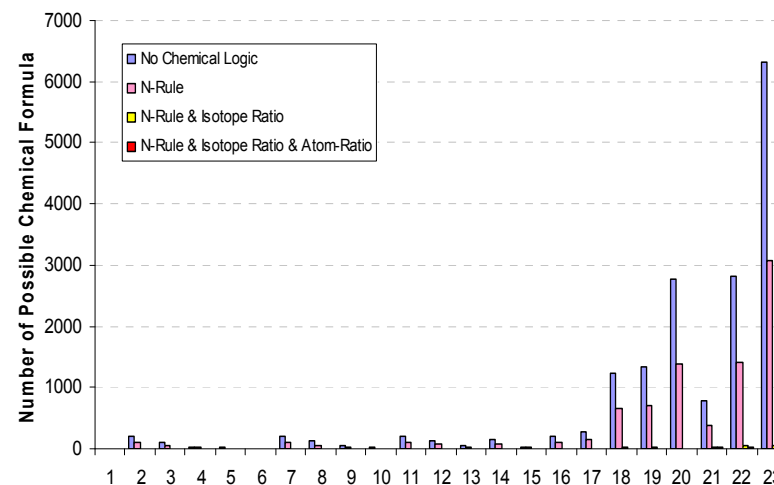

- Number of possible chemical formulas within 5ppm when considering
 - No chemical logic
 - N-rule
 - N-rule and Isotope ratio
 - N-rule, isotope and atom ratio

(considering limit of; C₅₀₀H₁₀₀₀O₁₀₀N₁₀₀P₅S₁₀F₁Cl₂Br₁)

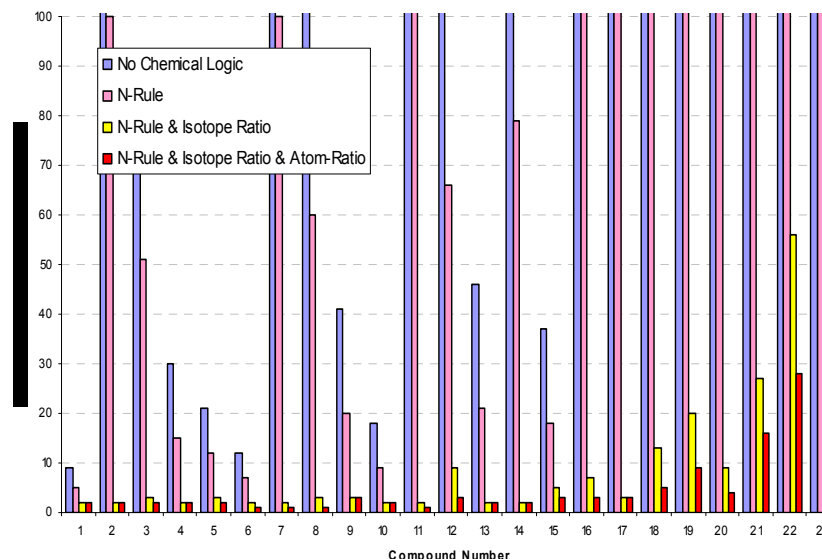




Identifying Biomarkers

- You have now found groups within the data, determined which masses are responsible for the groupings, and have elemental formulae for these masses
- Now what?
- Need to identify these compounds – many different metabolites can have the same elemental formula – you must determine structure for biomarkers
- Typical place to start is KEGG database, search for mass and elemental formula to try to identify compounds




How Much Mass Accuracy?

- Majority of compounds in database have molecular weights that are at least 10 ppm different from any other compound in database
- 3561 compounds in KEGG database are structural isomers, no amount of MS resolution or mass accuracy can separate these

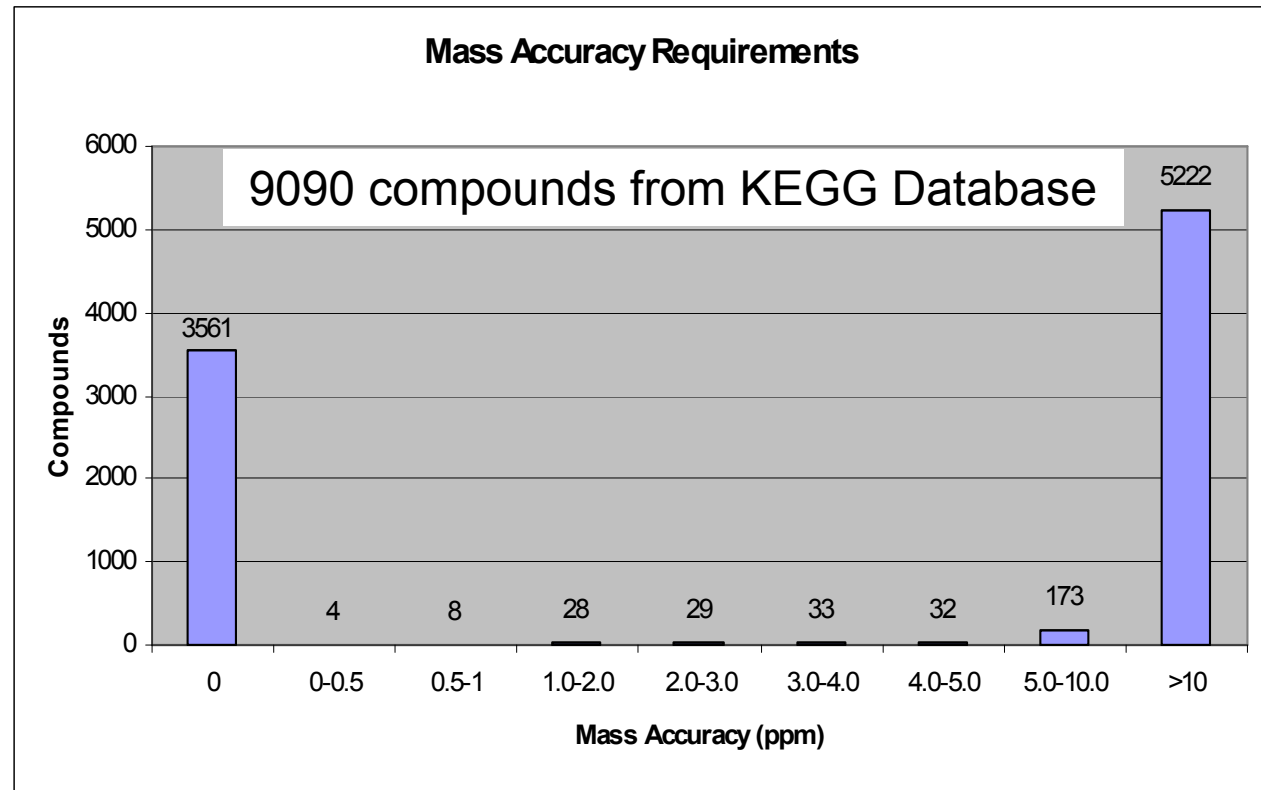

Sub-5 ppm mass accuracy gives only marginal improvement in compound identification when searching known databases

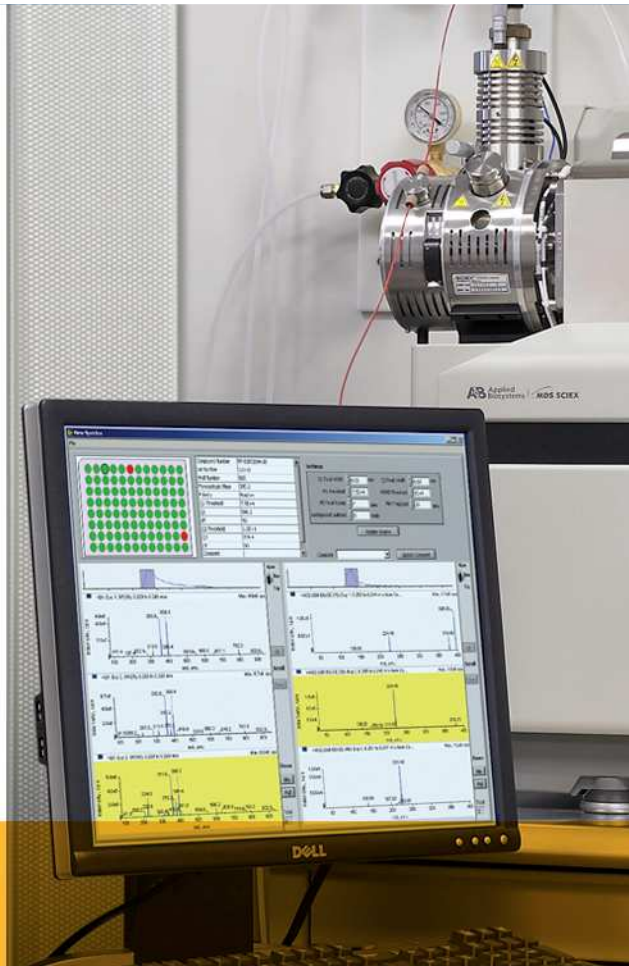

Targeted Metabolomics using Biocrates Absolute/IDQ™ kits





BIOCRATES' Approach To Metabolomics

Why Targeted Metabolomics?

- Robust and reproducible technology
- Absolute quantitation, not just relative peak intensities
 - Quantitative data is trusted data
 - (for you and regulatory authorities)
- Straightforward and rapid data interpretation because of identified and pre-annotated metabolites
- Closer to clinical needs due to resemblance to diagnostically established data formats (comparison to reference ranges)



AbsoluteIDQ™ Kit: Key Facts



- The AbsoluteIDQ Kit is designed to be used in combination with the API 4000™ or 4000 QTRAP® instrument
- Kit comes in 96-well plate format for both manual and automated operation
- Identifies and quantifies >150 different endogenous metabolites in blood plasma:
 - Acylcarnitines
 - Amino Acids
 - Glycerophospho- and sphingolipids
 - Hexose
- MS analysis by flow injection method, FIA-MS/MS (Quantification by MRM pairs)



Metabolite Details

163 Metabolites from four analyte classes

Acylcarnitines: 41

Amino acids: 13 proteinogenic + Ornithine

Phospho- and Sphingolipids

- Phosphatidylcholines 77 (Plasmalogene included)

- Lyso-Phosphatidylcholines 15

- Sphingomyelins 15

Hexose: sum of Hexose



MS instrumentation / robotic platform

Applied Biosystems:

Kit designed for Triple-quadrupole mass spectrometer from Applied Biosystems



API 4000™ or 4000 QTRAP®

Alliance with Hamilton Robotics:

Automated preparation possible (optional)
Help with integrating BIOCRATES methods on Hamilton robot will be provided




Hamilton Microlab Star




Kit Components

| Content | Description |
|---|---|
| Absolute/iDQ™ Kit Plate | Consists of a 96-well plate and filter plate attached with sealing tape |
| 96-Deep-Well Plate with Silicone Mat | For Autosampler Adjustment |
| BIOCRATES Solvent A | 250 ml MS running solvent |
| BIOCRATES Standards (Lyophilized) | 2 vials |
| BIOCRATES Quality Controls (QC) | 3 vials |
| Kit Manual | |
| USB Memory Stick | Has MetIQ Software and Oracle Express Database loaded. |


Unique Kit-Plate

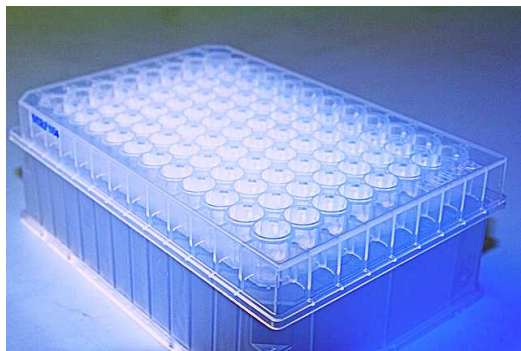

Proprietary Met/iQ™ Software

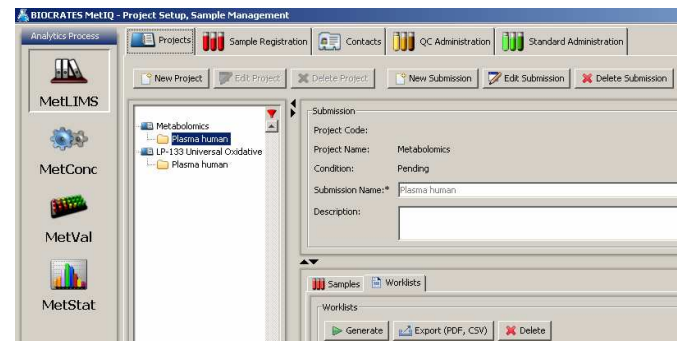

Overview of Workflow

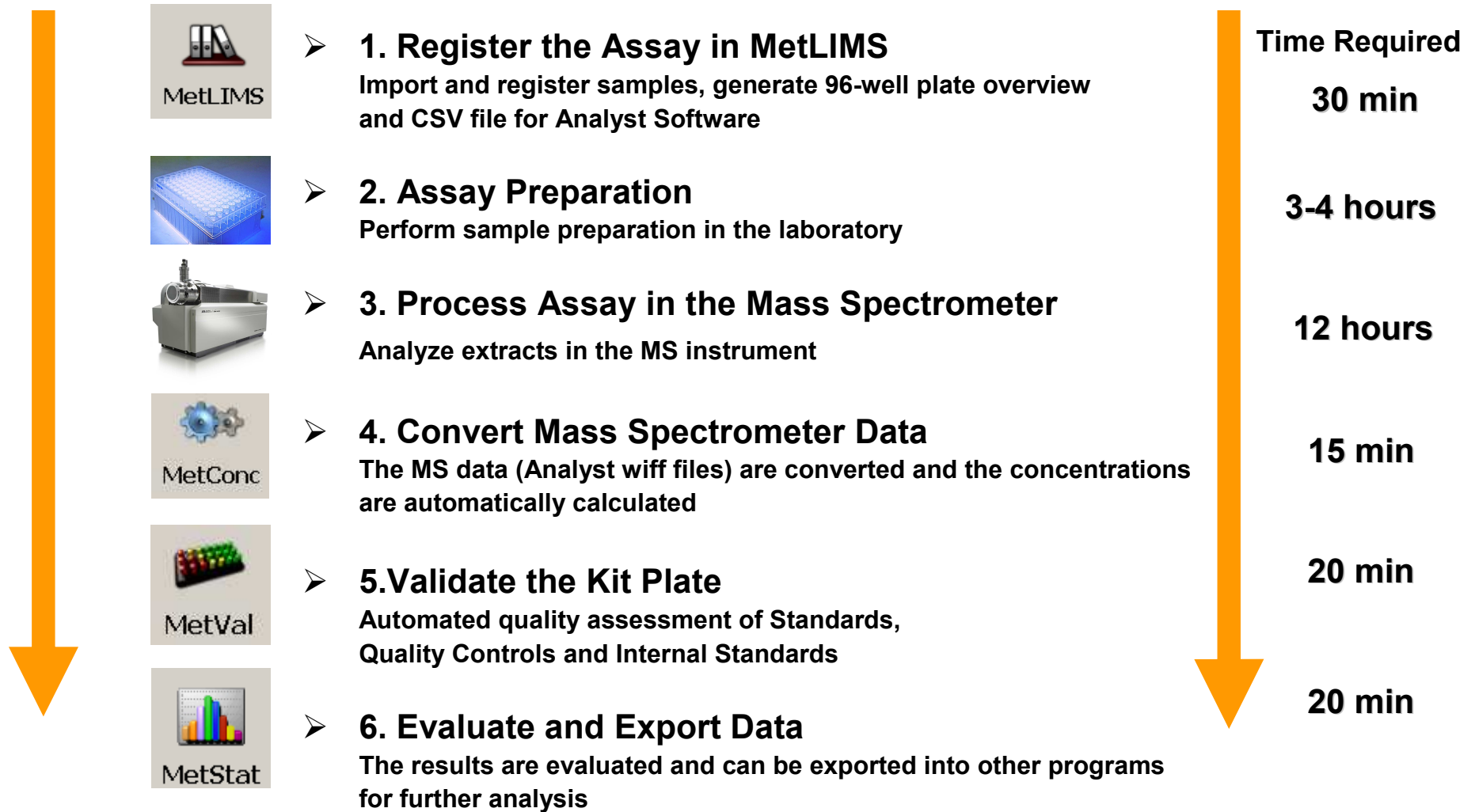




1. MetLIMS: Sample registration

96-well Plate Report and CSV File For Analyst

Two standards, three quality controls, one blank, one zero sample and the registered plasma samples are positioned on the kit plate

A 96-well plate report is printed out to help with the pipetting steps in the lab

A csv file is automatically generated for import into the Analyst acquisition batch

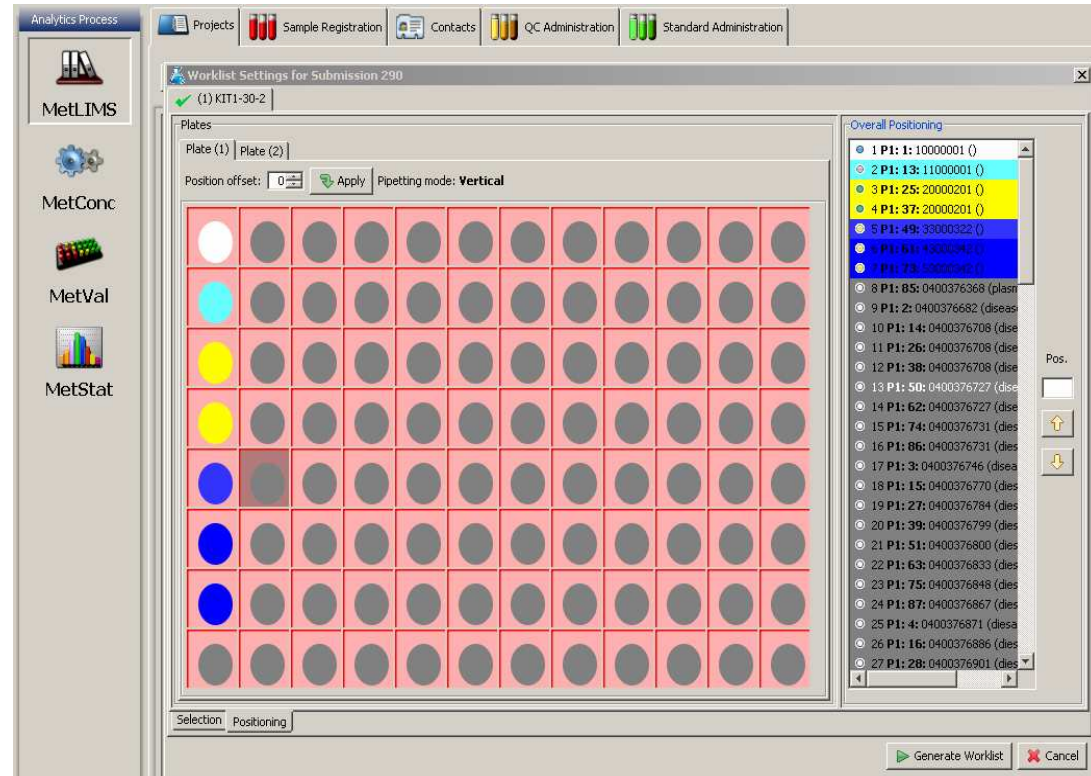




2. Assay Preparation: Unique Design of Kit Plate

➤ Advantages for sample preparation

**Upper Plate
For Extraction**

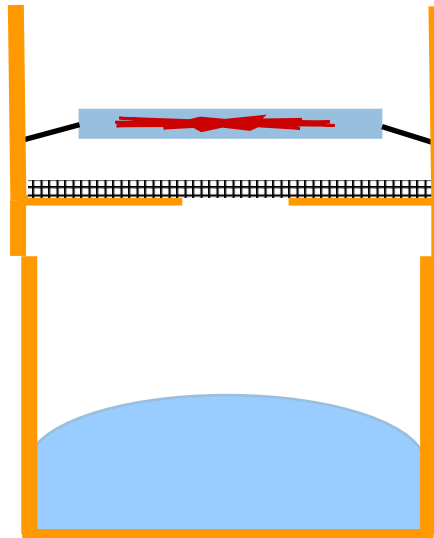


**Capture Plate
For Filtrate**

- **Internal standards incorporated**
- Protein precipitation and metabolite derivatization on upper filter
- Efficient circulation of nitrogen for drying step
- Efficient solvent extraction step



Very few sample preparations steps:

- 1) add 10 μL plasma (dry with nitrogen)
- 2) derivatize with PITC (dry with nitrogen)
- 3) extract metabolites (300 μL methanol)
- 4) dilute with MS running solvent

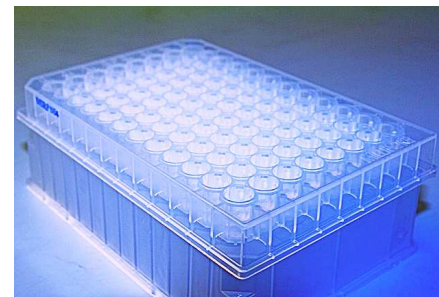

3. Process Assay in the Mass Spectrometer

175 MRM are measured in positive mode (22 IS)
 15 MRM are measured in negative mode (7 IS)
 2x 20 μ l injection, 30 μ l/min flow rate
 3 min run time, 7 min total time per sample

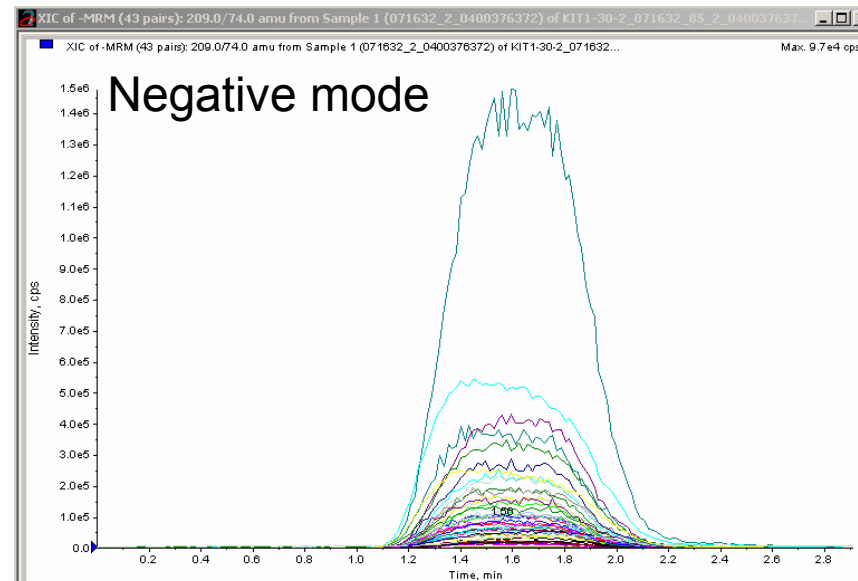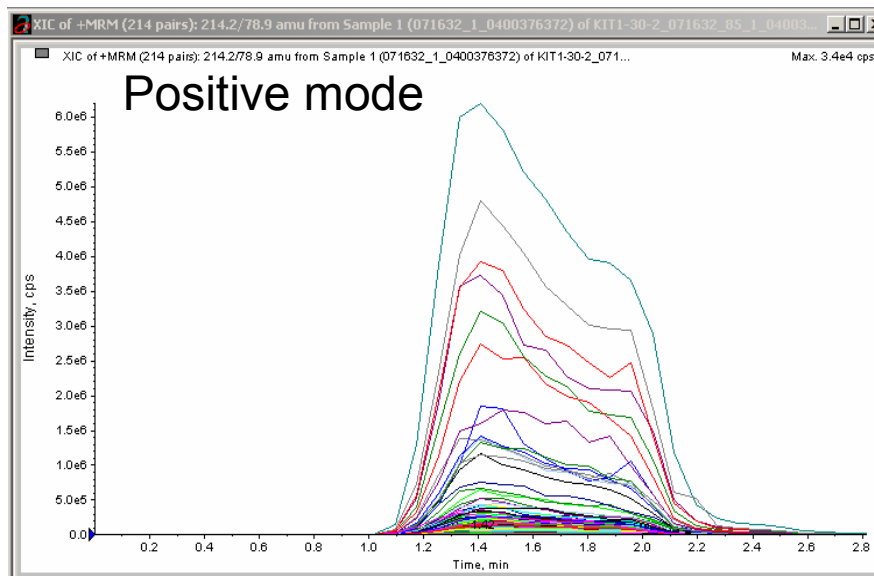


Calculation of concentrations is automated by Met/Q software




4. Convert Mass Spectrometric Data

- In the MetConc module the Analyst wiff files are converted and imported into the Met/Q Databank
- The concentrations of the metabolites are automatically calculated
- Direct link to Analyst software to open wiff files from this module

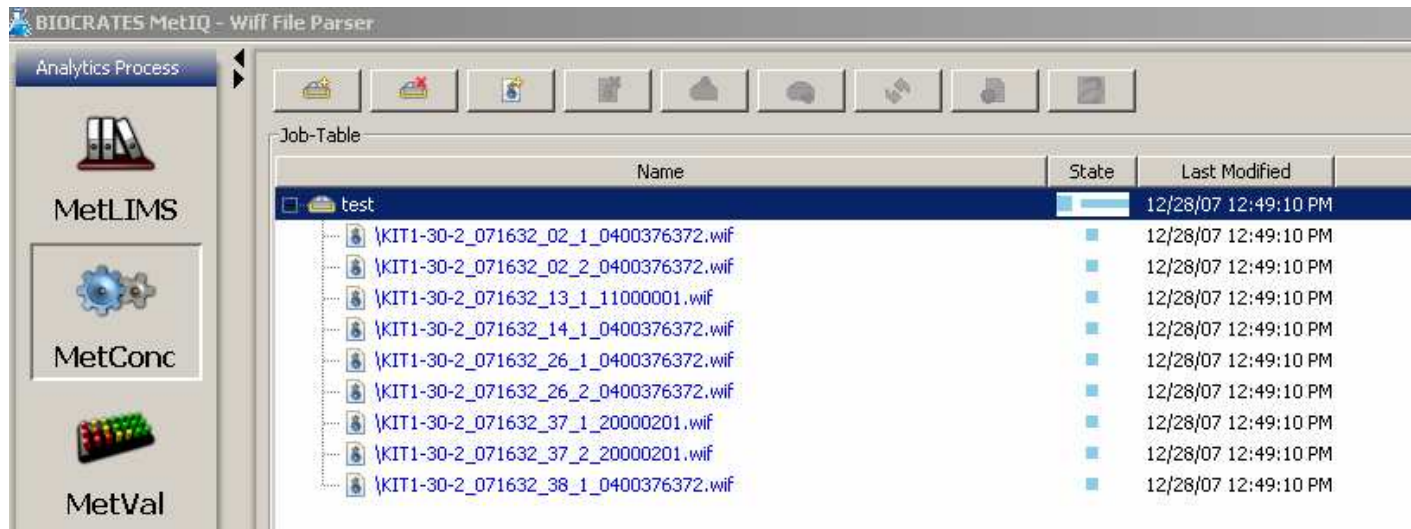

5. Validate the Kit Plate

- In the MetVal module an automated quality assessment of the Kit is performed
- Specifically, it is checked if the internal standards, standards, blank and quality control samples are within the specified ranges
- In case a value is out of range this will be indicated in red

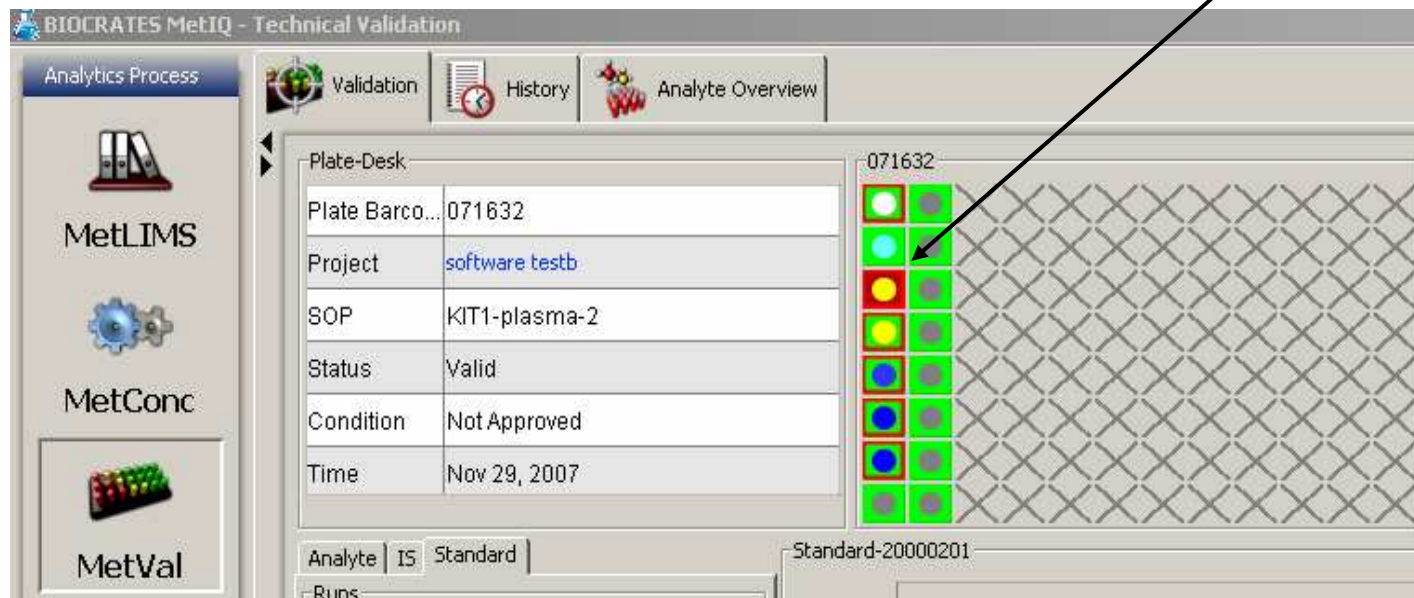

BIOCRATES MetIQ - Technical Validation

Analytics Process: **MetLIMS**, **MetConc**, **MetVal**

Validation | History | Analyte Overview

| Plate-Desk | |
|----------------|----------------|
| Plate Barco... | 071632 |
| Project | software testb |
| SOP | KIT1-plasma-2 |
| Status | Valid |
| Condition | Not Approved |
| Time | Nov 29, 2007 |


Analyte: IS Standard | Standard-20000201

Runs:



5. Graphical display of data in MetVal

Different graphs for easy visualization of validation results:
Standard graph: The BIOCRATES Standards (ES) are a mixture of analytes in identical concentrations as the Internal Standards (IS)
The graph shows the ratio of IS to ES – it should be close to 1

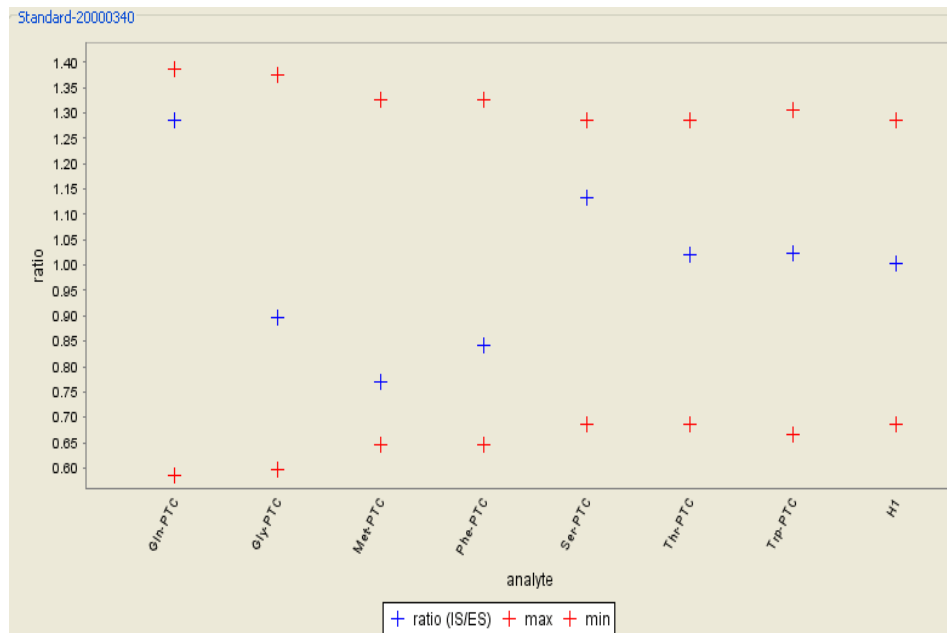

Red upper and lower range
Blue : measured values

6. Evaluate and Export Results

- The data are presented in tables sorted by options (analyte class, concentration, intensities)
- The information from the validation is transferred into the tables: In this example the blank value for the amino acid Thr was out of range as indicated by the red color
- The data can be exported from MetStat as a .txt or .csv file for import into other programs

BIOCRATES MetIQ - Data Extraction

Analytics Process | Link Samples | Display Data

MetLIMS | MetConc | MetVal | MetStat

| Met-PTC | Orn-PTC | Phe-PTC | Pro-PTC | Ser-PTC | Thr-PTC | Trp-PTC |
|------------|------------|------------|------------|------------|------------|------------|
| aminoacids | aminoacids | aminoacids | aminoacids | aminoacids | aminoacids | aminoacids |
| 9.067 | 21.289 | 74.258 | 1.935 | 2.292 | 7.007 | 2.48 |
| 222.290 | 79.335 | 97.209 | 218.452 | 152.297 | 189.223 | 74.11 |
| 18.280 | 3.147 | 4.425 | 0.656 | 2.309 | 9.855 | 1.05 |
| 139.754 | 64.749 | 75.228 | 202.565 | 124.695 | 159.536 | 63.94 |
| 263.231 | 68.984 | 73.607 | 56.228 | 70.094 | 62.433 | 53.63 |
| 116.264 | 64.639 | 75.452 | 202.286 | 121.703 | 149.607 | 63.72 |
| 219.647 | 65.143 | 73.517 | 56.309 | 70.811 | 62.185 | 52.21 |
| 104.975 | 56.845 | 73.793 | 197.613 | 118.227 | 146.504 | 61.80 |
| 150.727 | 63.115 | 75.106 | 195.829 | 120.905 | 160.311 | 62.38 |
| 105.377 | 65.029 | 75.885 | 201.585 | 123.455 | 150.451 | 63.67 |
| 148.704 | 59.866 | 73.745 | 191.487 | 115.466 | 151.461 | 60.66 |
| 106.304 | 59.577 | 76.438 | 205.303 | 121.594 | 154.485 | 64.06 |
| 162.822 | 60.733 | 72.499 | 196.799 | 117.668 | 152.371 | 60.51 |
| 109.292 | 61.840 | 73.789 | 197.632 | 120.040 | 157.878 | 62.37 |
| 182.060 | 57.872 | 69.638 | 186.217 | 113.120 | 139.875 | 59.81 |
| 150.064 | 61.285 | 76.861 | 204.273 | 122.159 | 166.492 | 64.18 |

Materials: plasma

Select the type of value to display:

- Concentration [µM]
- Intensity [Cps]
- Concentration Internal Std. [µM]
- Intensity Internal Std. [Cps]


Show class: aminoacids

Hide run numbers

Transpose values | Export



Variability in different laboratories

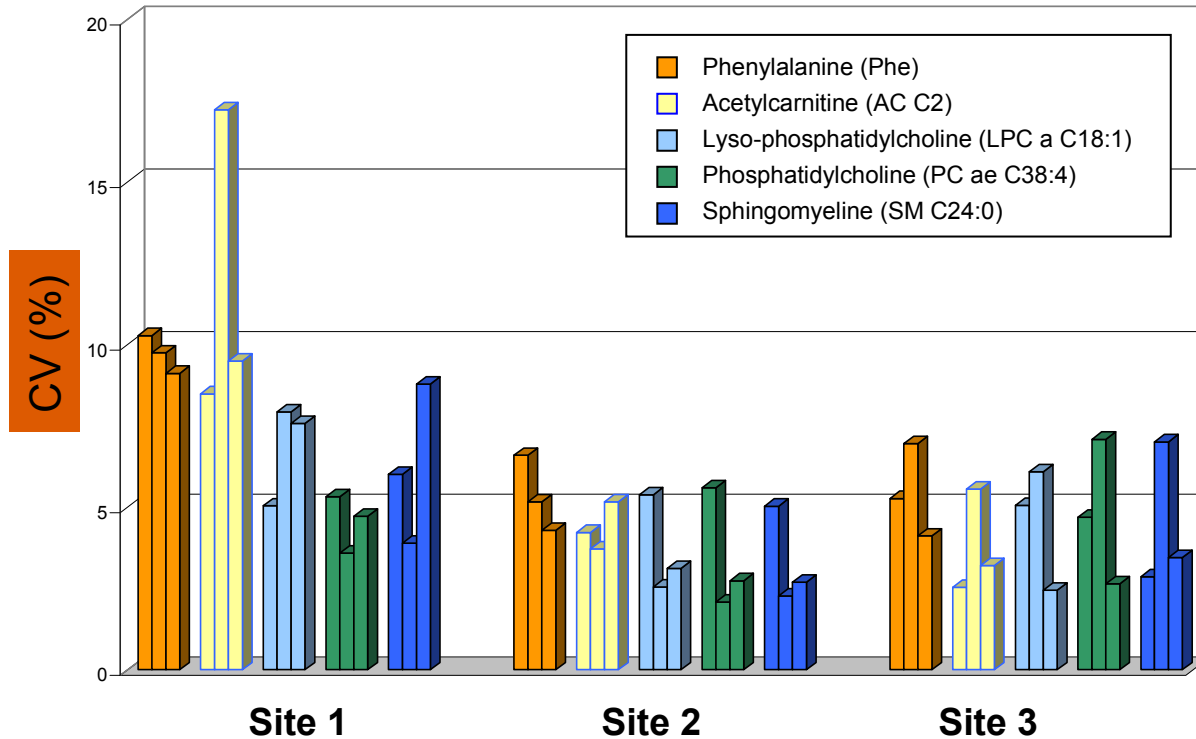

Representative results of test in three different laboratories:
 Coefficient of Variation (CV) of 5 metabolites at 3 different days and in 3 different sites are shown. In general the CVs were significantly below 15 % and comparable in the different labs.



Main Application Areas

Biomarker Research & Diagnostics

- Disease Phenotyping
- Metabolic Biomarker Discovery
- Early Disease Diagnosis


Pharmaceutical R&D

-Clinical & Pre-clinical studies

- Drug metabolism / Pharmacokinetics
- Pharmacodynamics / Efficacy
- Toxicity / Safety
- Quantitative comparison of drug candidates

Nutritional & Functional Food Analysis

- Health effects of functional food
- Effects of environment & exercise

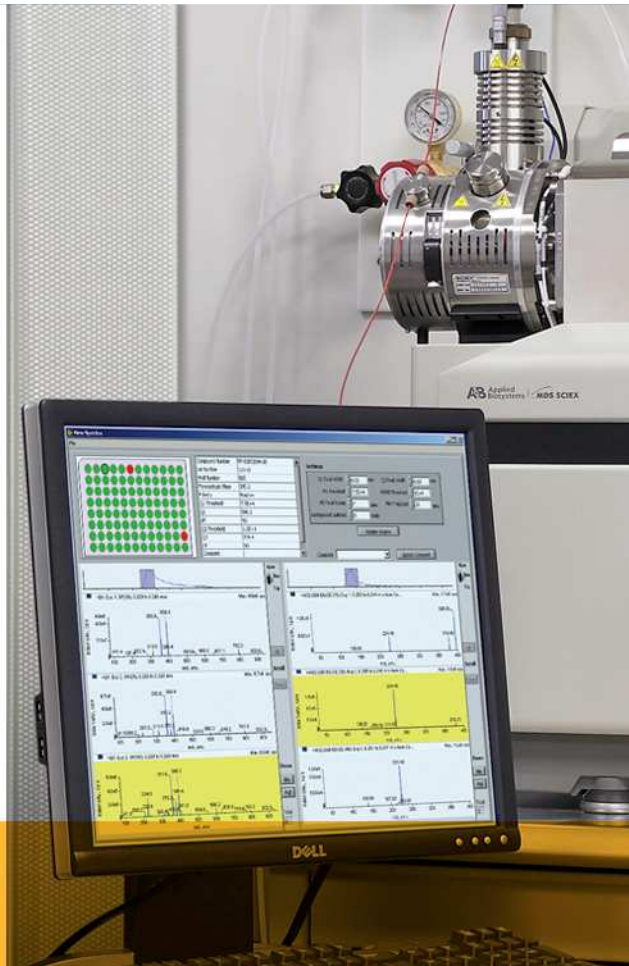


Lipidomics: *Applications & Workflows*



Lipidomics Overview: the “omics” era

Where does Lipidomics fit into the “omics” era?

- **Genomics & Proteomics** – mapping of genes, gene products and their interaction networks in a functioning cell
- **Metabolomics** – mapping of endogenous metabolites and how they’re affected by environmental conditions, including disease & toxicity, etc...
 - Metabolome is enormous and complex
 - Can only be approached in sections, e.g Lipids
- **Lipidomics** – characterization of the global changes in lipid metabolites (“lipidomics”)

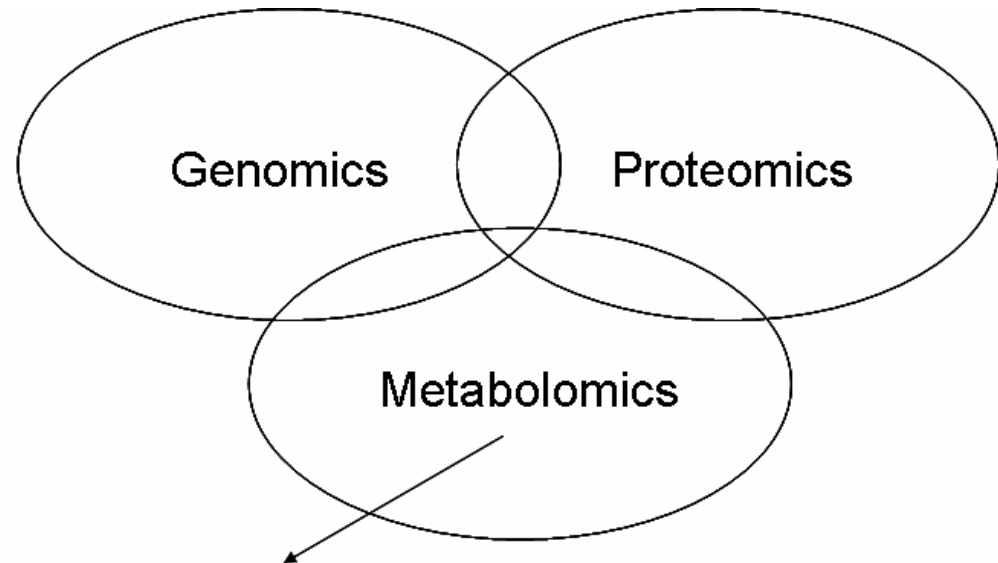


Approach in sections: eg. Lipids



Lipidomics Overview: What are they?

- Lipids are subdivided into classes:
 - Fatty Acids
 - Glycerolipids
 - Glycerophospholipids
 - Sphingolipids
 - Sterol Lipids
 - Prenol Lipids
 - Saccharolipids
- Nomenclature based on lipid class, head group and fatty acid chain length (with designations for number of double bonds – and location is potentially very important also)

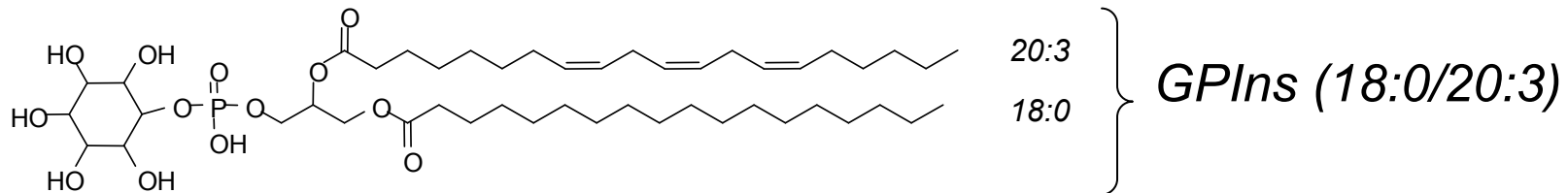

Glycerophospholipid inositol (GPIs)



Lipidomics Overview: Application Areas?

- **Pharma**
 - Better drug design (e.g. COX inhibitors)
 - Better drug delivery
 - Focus on obesity
- **Nutritional science**
 - Better health markers
 - Omega-3 fatty acids
- **Clinical – huge implications in a wide range of diseases**
 - Heart disease
 - Stroke
 - Cancer
 - Alzheimer's
 - Possible to have a diagnostic tool around specific lipid monitoring (lipid biomarkers)




Software Enabling Lipidomics Workflows

LipidProfiler™ (*Prototype Research Grade Software*)

- *Lipid Profiler* is a data processing tool which identifies and semi-quantitates lipids from electrospray MS and MS/MS data.
- Three modes of data processing:
 1. information for all m/z ion peaks
 2. m/z ion peaks for a set of target lipid species
 3. m/z masses for characteristic lipid masses (Identify species mode)
- Qualitative profiling of lipid MS data is validated by links to the information in the database for characteristic lipid headgroup, fatty acid, or long chain base fragments and neutral losses.
- The peak profiles, lipid profiles, and fatty acid profiles are displayed in a graphical interface which can be exported from the application along with all numerical quantitative data.
- Compatible with Analyst® QS and Analyst® Software



MarkerView™ and LipidProfiler™ Workflow:

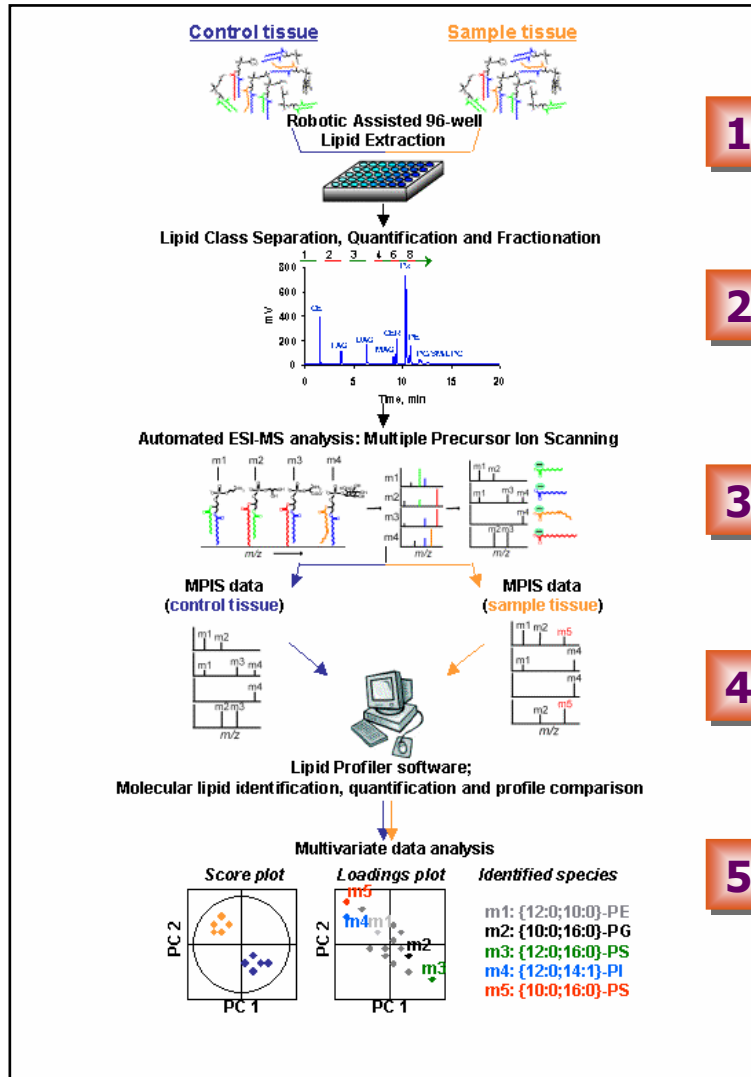

1. Robotic Assisted 96-well Total Lipid Extraction
2. Lipid Class Separation, Quantification & Fractionation
3. Robotic Sampling and Automated MS Acquisition
4. Computer-assisted Identification & Quantification of Molecular Lipids (*LipidProfiler*)
5. Multivariate & Univariate Data Analysis (*MarkerView*)

Kim Ekroos, VTT, Finland

Workflow

Load MS Data

Processing Setup

Find Lipids

Results

Preview Report

Summary

4 files, 10 samples

4 methods used

Results Available

0 snapshots

5 pages in report

LP Utilities

Catalogue

Calculator

Results

Go to Next

View Mode:

Peak Intensities

Grid...

Column: File 'Lipid2.wiff', Sample 'SampleName1', Period 1. Path = C:\Analyst\Projects\AnalystProject1\Lipid2.wiff

Options:

Sort by Lipid ID

Sort by Count

Show Experiment Details

Cumulative Selections

Export Data...

Filter Data:

All

Profile Tests...

Samples...

| Lipids \ Samples-Periods | Count | Smp1, Prd1 | Smp1, Prd1 | Smp2, Prd1 | Smp2, Prd2 | Smp1, Prd1 | Smp2, Prd1 | Smp2, Prd2 | Smp3, Prd1 |
|--------------------------|-------|------------|------------|------------|------------|------------|------------|------------|------------|
| DAG 31:4+NH4 (DAG 14:1) | 14 | 274.177 | 156.148 | 0.000 | 22.785 | 410.213 | 37.951 | 433.508 | 196.1 |
| DAG 32:1+NH4 (DAG 14:0) | 12 | 125.803 | 490.322 | 0.000 | 67.483 | 343.269 | 0.000 | 64.425 | 0.0 |
| DAG 31:4+NH4 (DAG 16:2) | 13 | 0.000 | 55.066 | 454.373 | 0.000 | 199.389 | 0.000 | 321.763 | 223.0 |
| DAG 34:3+NH4 (DAG 16:2) | 12 | 192.658 | 91.341 | 140.518 | 278.362 | 0.000 | 332.778 | 204.988 | 0.0 |
| DAG 32:1+NH4 (DAG 16:1) | 12 | 235.837 | 88.591 | 465.532 | 59.386 | 210.492 | 254.208 | 0.000 | 0.0 |
| DAG 34:3+NH4 (DAG 16:1) | 18 | 139.227 | 37.748 | 115.712 | 300.896 | 465.095 | 278.784 | 0.000 | 214.0 |
| DAG 34:2+NH4 (DAG 16:1) | 14 | 352.516 | 0.000 | 278.898 | 277.802 | 262.389 | 0.000 | 253.078 | 445.0 |
| DAG 32:1+NH4 (DAG 16:0) | 13 | 0.000 | 199.729 | 73.859 | 0.000 | 0.000 | 0.000 | 273.983 | 120.0 |
| DAG 34:2+NH4 (DAG 16:0) | 14 | 74.830 | 184.351 | 0.000 | 0.000 | 0.000 | 480.287 | 255.278 | 121.0 |

Profiles

Data: Details

Data: Spectra

Add to Report

Main Plot Type

Peak Profile

Show 2nd Plot

2nd Plot Type

Molecular Species

Plot Attributes

Show Legend


Logarithmic Y Scale

Group Sample Data

Group Lipid Data

Normalize for Sample

Show Statistics in Grid

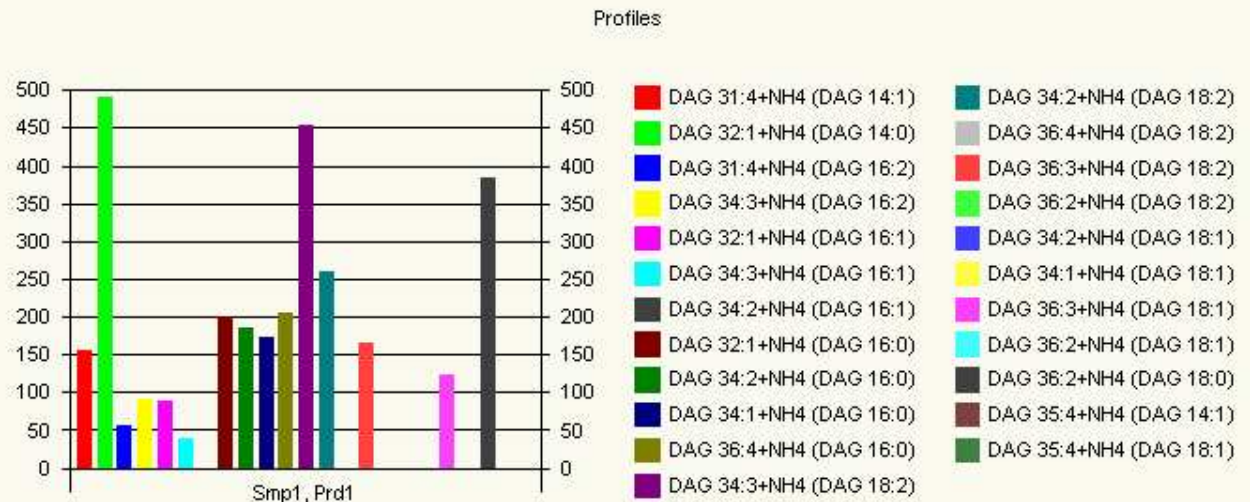

| Selection | Smp1, Prd1 |
|-------------------------|------------|
| DAG 31:4+NH4 (DAG 14:1) | 156.148 |
| DAG 32:1+NH4 (DAG 14:0) | 490.322 |



Lipid Species Profile



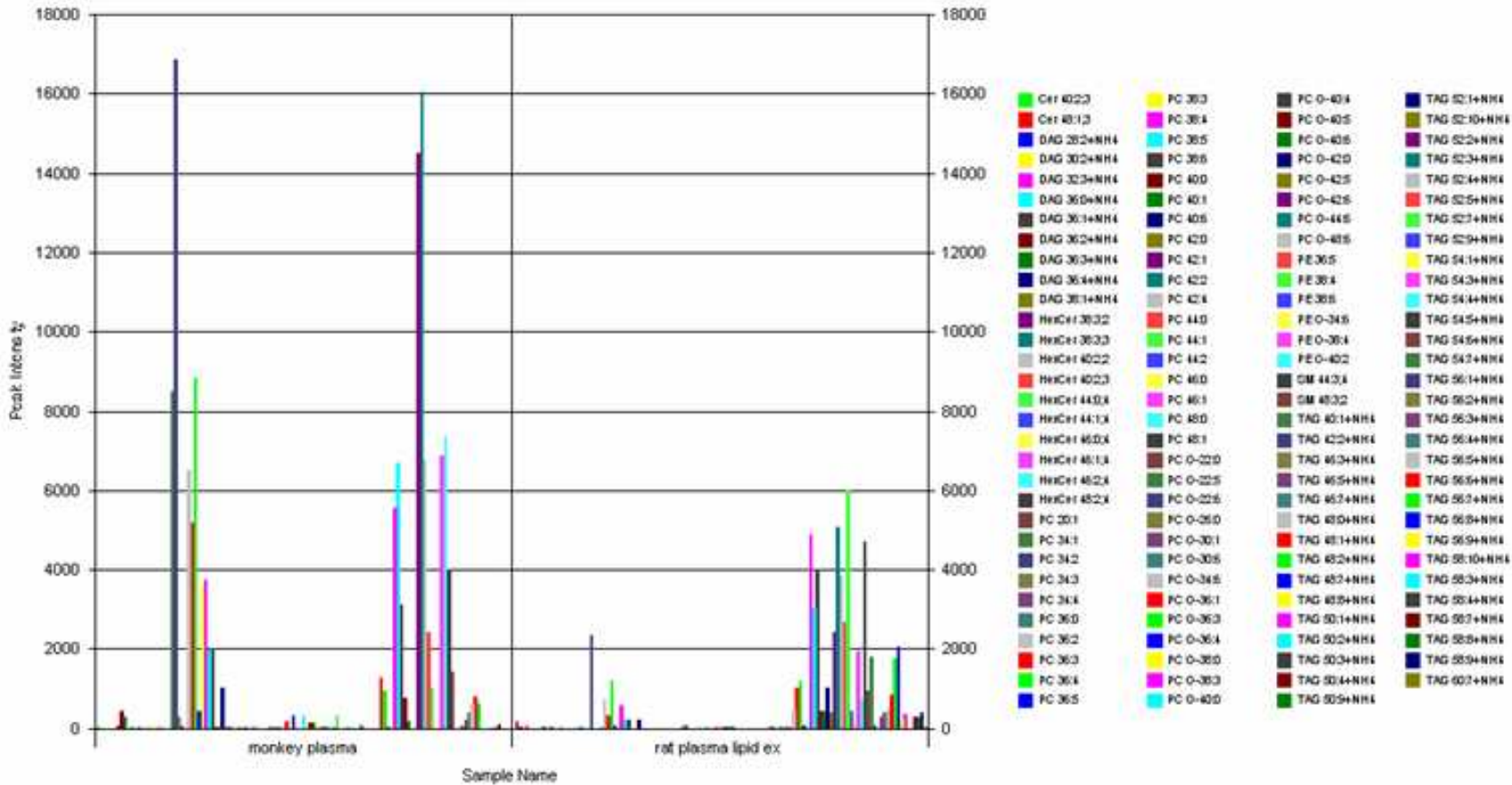


Figure 3. Lipid species profile comparisons of monkey and rat plasma lipid extracts generated by processing QSTAR[®] Elite +TOF MS and IDA product ion data through LIPID PROFILER. Over 120 lipid species were identified through a glycerophospholipid database targeted analysis.

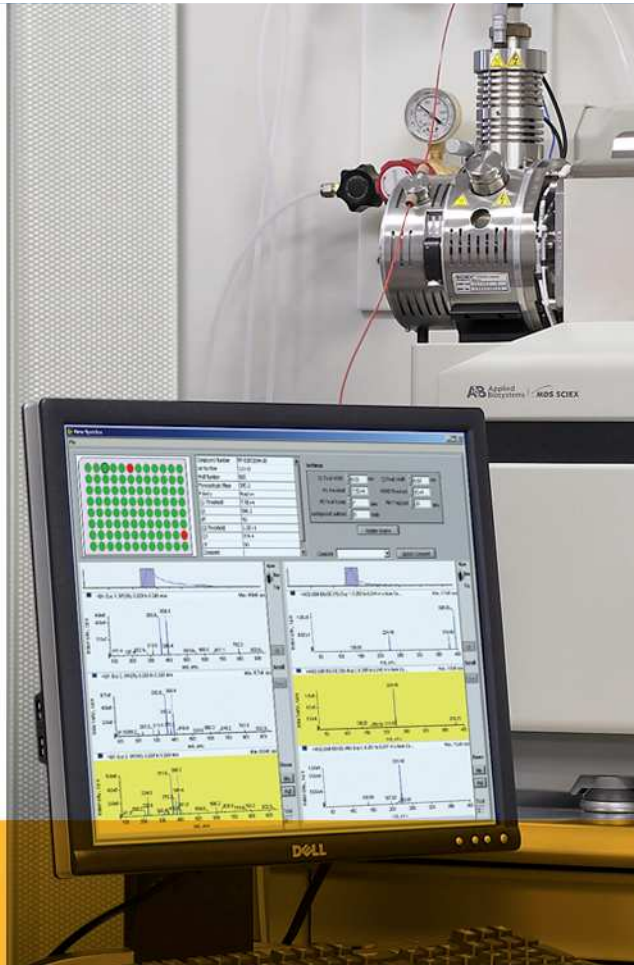

Thank You!
Questions?

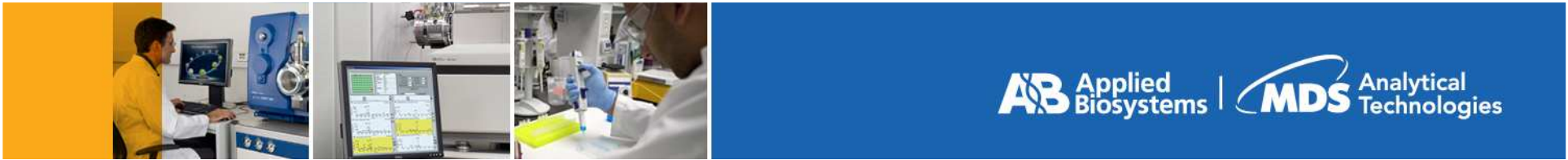


Trademarks/Licensing

For research use only. Not for use in diagnostic procedures.

AB (Design) and Applied Biosystems and are registered trademarks of Applied Biosystems Inc. or its subsidiaries in the US and/or certain other countries.

MIDAS, API 5000, AcQuRate, eQ, DiscoveryQuant, iMethods, MRMPilot, ProteinPilot, MultiQuant and AB SCIEX Triple Quad are trademarks and QTRAP, Cliquid, LightSight, and LINAC are registered trademark of Applied Biosystems/MDS Analytical Technologies, a joint venture between Applied Biosystems Inc and MDS Inc.

© 2008 Applied Biosystems Inc. and MDS Inc. Joint Owners. All rights reserved.

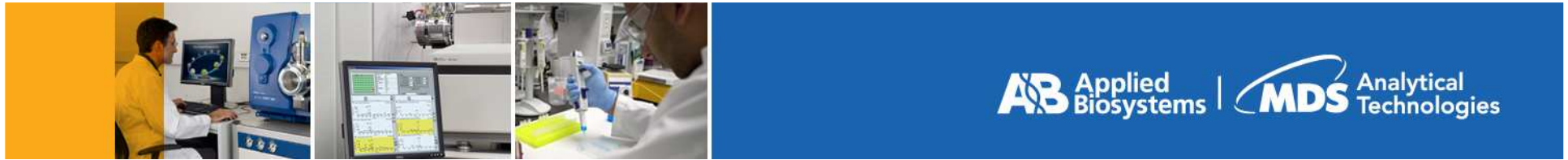

For Research Use Only. Not for use in diagnostic procedures.

Applera, Applied Biosystems and AB (design) are registered trademarks of Applera Corporation or its subsidiaries in the US and/or certain other countries.

MarkerView is a trademark and Analyst, QTRAP and QSTAR are registered trademarks of Applied Biosystems/MDS Analytical Technologies, a joint venture between Applera Corporation and MDS Inc.

All other trademarks are the sole property of their respective owners.

© 2008 Applera Corporation and MDS Inc. Joint Owners. All Rights Reserved.